

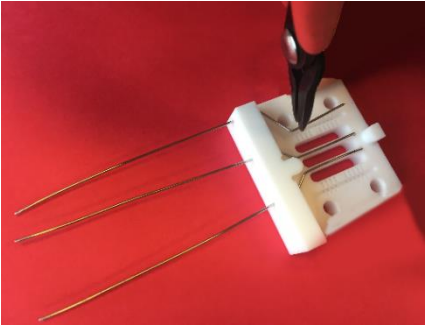

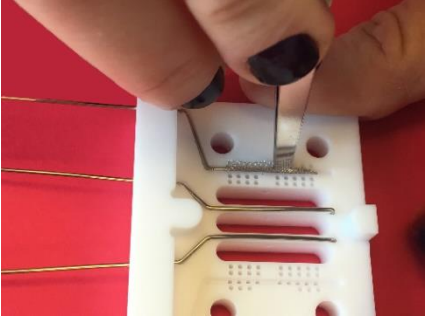


Instructions: Pt mesh replacement in a BekkTech BT-11x (-110, -112, -115, or -710)

** Note: Pre-cut replacement mesh is available from Scribner - contact us for a quote **

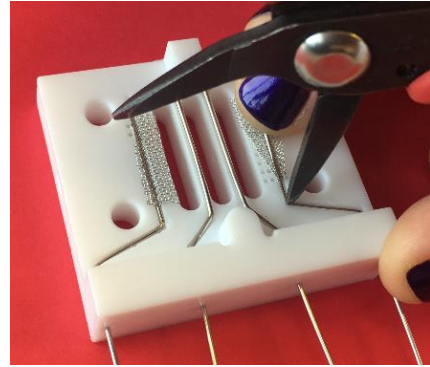
Required Parts:

1. Replacement Pt mesh
2. Tweezers or needle-nose pliers
3. Small flat-head screw driver

| | |
|--|--|
| <p><i>CAREFULLY</i> pry the Pt wire from the groove with a small, flathead screwdriver, needle-nose pliers or tweezers</p> <p>NOTE: There is no need to remove the wire entirely from the clamp; the wire only needs to be popped out from the groove</p> |  |
| <p>Slip the old mesh off the Pt wire</p> <p>Use tweezers to hold the replacement Pt mesh piece over the grooved clamp (area with the tiny holes) and under the Pt wire</p> |  |
| <p>Push the wire and mesh into the groove using a small screwdriver or tweezers</p> |  |



Use needle-nose pliers to firmly press the wire into the groove in several locations simultaneously



Carefully flatten the Pt mesh against the clamp using tweezers

